


Navigator Login Screen

- Use your provided login and password to access Navigator.
- If you forget your password, click the **Forgot Password** option below **LOGIN**

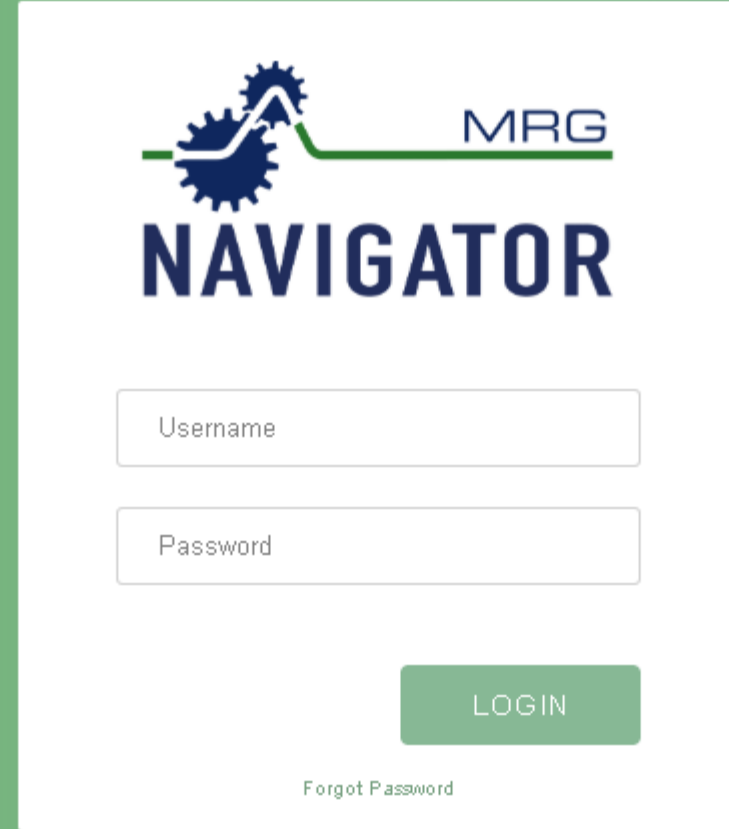

MRG
NAVIGATOR

Username

Password

LOGIN

[Forgot Password](#)

Welcome! To view your sample results, enter your User ID and Password then click the Login button. If you need assistance or additional sampling supplies, please contact your MRG representative, or call the lab at 717-843-8884. To begin with new sample information, click the New button under the Barcoded tab.

Navigator Main Screen

1. The default view of samples will be the “Grease” dashboard
2. Hover over “SAMPLES” to switch screens to “OIL”
3. Use accounts to view account information and change Password

The screenshot displays the Navigator Main Screen interface. At the top left, the logo for 'MRC NAVIGATOR' is visible. A navigation menu on the left side includes 'SAMPLES' (highlighted with a blue hexagon labeled '2') and 'WEBVIEW'. The main content area is titled 'Grease Samples' (indicated by a blue hexagon labeled '1') and features four summary cards: 'Recorded' (0), 'Being Analyzed' (0), 'Ready For You' (0), and 'History' (0). A 'NEW' button is located below the cards. A search bar with a 'FILTER' dropdown and a search icon is positioned to the right. Below the search bar, a table header lists various columns: SAMPLE ID, COMPANY, EQUIPMENT TYPE, EQUIPMENT ID, SAMPLE POINT, IDENTIFIED BASELINE, RUNTIME, FACILITY, AREA, ROUTE, APPLICATION, and INTEI. The table content area is currently empty, displaying the message 'No data available in table'. In the top right corner, there is a 'Log Out' button and an 'Account' link (highlighted with a blue hexagon labeled '3').

Accounts

1. In addition to changing your password, you can update how you would like to receive information from the system.
2. Set email reminders when you have to sample machinery and also set the interval date frequency
3. Have reports emailed to you. Reports will send as one pdf document versus individual documents of samples in **Ready For You** status.

Edit Notifications

Email Reminders

Receive email reminders to resample your machinery

Reminder Interval

How many days after you complete your samples until you receive a reminder

Email Reports

Reports will be emailed to you after analysis is complete

CANCEL

SAVE

Recorded Screen

1. Click on new and follow the New Sample Input Directions to enter sample information. Several drop-down options are available to make sure samples are entered in consistently.
2. Samples entered will show up in this area. Once entered, fields may be edited from table format or the edit screen

The screenshot displays the MRG NAVIGATOR interface for Oil Samples. The top navigation bar includes the MRG NAVIGATOR logo, a back arrow, the title 'Oil Samples', and links for 'Account' and 'Log Out'. The left sidebar contains 'SAMPLES' and 'WEBVIEW' options. The main content area features four summary cards: 'Recorded' (0), 'Being Analyzed' (0), 'Ready For You' (3), and 'History' (0). A 'NEW' button is located below the cards, highlighted with a blue circle and the number 1. Below the cards is a table with columns for SAMPLE ID, COMPANY, EQUIPMENT TYPE, EQUIPMENT ID, SAMPLE POINT, IDENTIFIED BASELINE, RUNTIME, FACILITY, AREA, ROUTE, APPLICATION, and INTERVAL. The table is currently empty, showing 'No data available in table' and 'Showing 0 to 0 of 0 entries'. A blue circle with the number 2 is positioned over the table area.

Being Analyzed Screen

1. Once samples are received, they move to **Being Analyzed**
2. Use **EDIT** to update sample information if anything changes from initial entry
3. Use the **FILTER** options to show/hide entry and sample fields

The screenshot displays the 'Oil Samples' dashboard. At the top left is the title 'Oil Samples', and at the top right are links for 'Account' and 'Log Out'. Below the title are four status cards: 'Recorded' (8 items, blue icon), 'Being Analyzed' (1 item, red icon), 'Ready For You' (6 items, green icon), and 'History' (0 items, yellow icon). Below these cards are five buttons: 'NEW', 'EDIT', 'BASELINE', 'IMAGES', and 'DELETE'. To the right of these buttons is a search bar with a magnifying glass icon and a 'FILTER' button. Below the buttons and search bar is a table with columns for 'SAMPLE ID', 'SAMPLE DATE', 'EQUIPMENT TYPE', 'EQUIPMENT ID', 'SAMPLE POINT', 'IDENTIFIED BASELINE', 'RUNTIME', 'FACILITY', 'AREA', 'ROUTE', 'APPLICATION', and 'INTERVAL'. The first row of data shows 'Oii57', 'Type', 'ID', 'Point', 'Baseline', '2000', 'Fac', and 'App'. Blue callout boxes with numbers 1, 2, and 3 are overlaid on the interface: '1' points to the table header, '2' points to the 'EDIT' button, and '3' points to the 'FILTER' button.

Oil Samples Account Log Out

8 Recorded 1 Being Analyzed 6 Ready For You 0 History

NEW EDIT BASELINE IMAGES DELETE

FILTER

SAMPLE ID	SAMPLE DATE	EQUIPMENT TYPE	EQUIPMENT ID	SAMPLE POINT	IDENTIFIED BASELINE	RUNTIME	FACILITY	AREA	ROUTE	APPLICATION	INTERVAL
Oii57		Type	ID	Point	Baseline	2000	Fac			App	

Ready for You Screen

1. View data in a table format. Like the **Being Analyzed** page, filters may be set to limit data fields observed.
2. Download an excel file of the data in the current view
3. Print individual or multiple PDF reports for samples
4. Once Data is reviewed, send the sample to **History** for archive purposes. This keeps **Ready for You** for samples which need reviewed

Oil Samples

Companies

8 Recorded

1 Being Analyzed

6 Ready For You

0 History

SELECT ALL EXCEL REPORT HISTORY BASELINE IMAGES EDIT DELETE


SAMPLE ID COMPANY EQUIPMENT TYPE EQUIPMENT ID SAMPLE POINT IDENTIFIED BASELINE CORRECTED BASELINE RUNTIME FACILITY AREA ROUTE APPLICATION


SAMPLE ID	COMPANY	EQUIPMENT TYPE	EQUIPMENT ID	SAMPLE POINT	IDENTIFIED BASELINE	CORRECTED BASELINE	RUNTIME	FACILITY	AREA	ROUTE	APPLICATION
Oil17	Demo	Motor	Motor23	Drain	Example Baseline	Example Baseline	900	Bld1	Area	Route	Hvac
Oil30	Demo	Type	ID	Point	Example Baseline	Example Baseline	2000	Facility			App
Oil60	Demo	Motor	Motor1	Drain	AW-46	AW-46	600	Bld1			Hvac
oil4	Demo				AW-46	Mobil SHC 630 2240.54	5				


History Screen


1. Samples sent to History will display here.
2. All the features for reports/excel functions are the same as in the **Ready For You** screen.

Oil Samples Account [Log Out](#)


8
Recorded


1
Being Analyzed


6
Ready For You


0
History

SELECT ALL EXCEL FILTER

SAMPLE ID	COMPANY	EQUIPMENT TYPE	EQUIPMENT ID	SAMPLE POINT	IDENTIFIED BASELINE	CORRECTED BASELINE	RUNTIME	FACILITY	AREA	ROUTE	APPLICATION
-----------	---------	----------------	--------------	--------------	---------------------	--------------------	---------	----------	------	-------	-------------